Modelo 146 P | [Model 146P](#)

Marca | [Trademark: Zuzi](#)

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Zuzi microscope 146 P is equipped with infinity corrected optics. It gives excellent technical features in samples observation under polarizing light. It is very useful in petrography, mineralogy, materials chemistry and in histology and pathological anatomy to observe the numerous structures that have birefringence (estructuras cristalinas, pigmentos, lípidos, depósitos de amiloide, etc.).

For that reason, the microscope has a polarized filter under the slide that transmit the polarized light in direction N-S, and a analyzer filter, settle on top of the slide, that transmit only the light that is vibrating in direction E-O. When both polarizes are in orthogonal position (placed with 90° between each other), the light generated by the polarizer is extinguished and you can not see anything; however, if the sample shows birefringence the birefringence structure is illuminated over a dark background.

Zuzi microscope 146P present infinity corrected optic and lighting system that guarantee to obtain high quality images free of distortions and mistakes.

The equipment has a Köhler lighting system with field diaphragms and a opening; graduated precondenser and Abbe condenser (AN:1.25) adjustable in height by a pinion and rack. The lighting is made by a halogen lamp of 30W with the regulator of intensity. The objectives, quality achromatic planes, are centred in the nosepiece, avoiding deviations in the light.

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The equipment has a polarized filter under the condenser that can be turn in a 360 ° angle and it is graduated. The analyzer filter, settle in the intermediate tube between the nosepiece and the head, presents a rotation angle of 90° and a Vernier graduation.

The intermediate tube has a Bertrand lens easily slider from the optic way to conosopic observations and a crack for the insertion of compensative plaque between the ocular and the objective. The equipment has some accessories wedges, a gypsum plaque (1 l), mica plaque (1/4 l) and a quartz wedge that are inserted at 45° of the vibration directions of the polarizer and the analyzer.

The stage, rotary and circular, show a 160mm diameter and a graduation from 0° to 360° with Vernier giving a resolution of 0.1° and a good accuracy in the determination of the angles rotation. A mechanism lets fix the stage in many positions.

The slide has as well two clamps to hold the sample and the possibility of joining a mechanic cart that gives a comfortable and quick moving of the sample and the location of the interested structures.

As important as technical features, is the ergonomic design that assures a natural position of the head and the body to avoid fatigue when the microscope is going to be use for a long time. For that reason, the microscope 146P has a Siedentopf binocular head inclined 30° with regulation of the interpupilar distance and the dioptrically correction, equipped with oculars suitable for users with glasses.

The controls macro and micro coaxial are in a low position to obtain a comfortable use. The right control have a control ring of friction and the left one has a block lever for the higher limit of focusing to avoid the samples hit accidentally the lens.

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